Application Note for LabVIEW Programming

Summary

1.	Pre	face	3
2.	Inst	ructions	3
	2.1.	Preliminary verifications	3
2	2.2.	Download VIs and unzip files	3
2	<u>2.3.</u>	Create a LabVIEW project	4
2	2.4.	Create your first LabVIEW program	6
3.	Erro	Drs	11
4.	Con	nmunication issues	11

1. Preface

This document will help you to start a project with LabVIEW in order to drive our product like:

- CCS/std
- CCSI
- Central / Central I
- PDM HPP
- Tombak

2. Instructions

2.1. Preliminary verifications

Please verify the correct functioning of the product before to start a LabVIEW project.

2.2. Download VIs and unzip files

The first step is to unzip the files and place them in this specific folder:

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Figure 1 : Unzipping the files in the correct folder

2.3. Create a LabVIEW project

1- Open LabVIEW and Select "create a project"

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2- Add the Open_Communication.vi and Open_Communication.vi, these are the first VIs to add in order to drive our product

Or you can directly add the VI to the project.

3- You can press CTRL-H to open a help window

Thanks to this window we can know the input ant output signals. In this Open_Communication.vi, there are 2 input signals (portName / input error) and 3 output signals (portPointer / function return / output signal).

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5- Check the portName (USB serial Port) with the device manager.

Ports (COM et LPT)
USB Serial Port (COM41)

In this case, the portName is COM41. So, we must write "COM41" in the input signal portName.

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If there is no problem, the "function return" signal returns 0. The portPointer is generated by the Open_Communication.vi

7- Add different VIs from our library to drive the product

This is an example of using our LabVIEW function, we just set and read the temperature. **Don't forget** to use the Apply_Request.vi after Set_XXX.vi (for example: set_temperature.vi or set_pulse_width.vi) and the Close_Communication.vi at the end of your program !

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3. Errors

RS232 error codes :

- 0 (SP_OK) : Operation completed successfully.
- -1 (SP_ERR_ARG) : Invalid arguments were passed to the function.
- -2 (SP_ERR_FAIL) : A system error occurred while executing the operation.
- -3 (SP_ERR_MEM) : A memory allocation failed while executing the operation.
- -4 (SP_ERR_SUPP) : The requested operation is not supported by this system or device.

Laser error codes :

0 : No error -30001 : Timeout Error -30002 : Unknown Command -30004 : Request Error -30008 : Length error -30016 : Checksum Error

4. Communication issues

If communication problems appear check the following points :

1) With the device manager check that you enter the correct Com Port

- 2) Please check if the GUI is not connected to the product
- 3) Be sure to put the close_com_vi at the end of your program. Otherwise LabView will keep the Port COM open and the new process will not be able to connect to the same Port Com. In this

case, disconnect and reconnect the USB/JACK cable. If the problem persists, remember to check that the drivers are up to date on the official FTDI website: https://ftdichip.com/drivers/vcp-drivers/

4) If the program crashes during the execution, the Port COM is still open. So disconnect and reconnect the USB/JACK cable